


200 MHz Variable Gain Photoreceiver

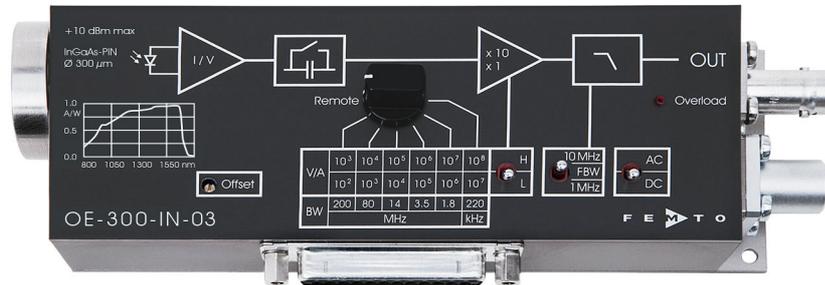

The image shows model OE-300-IN-03-FST with 1.035"-40 threaded flange and coupler ring.


Features

- **Adjustable transimpedance gain from 10^2 to 10^8 V/A**
- **Wide bandwidth up to 200 MHz**
- **InGaAs-PIN photodiode covering the 800 to 1700 nm wavelength range**
- **High dynamic input range up to 10 mW optical power**
- **Very low noise, NEP down to 52 fW/√Hz**
- **Switchable low pass filters for minimizing wideband noise**
- **Threaded 1.035"-40 and unthreaded 25 mm dia. free space input available, compatible with many optical standard accessories**
- **Full manual and remote control capability**

Applications

- **All-purpose low-noise photoreceiver (O/E converter) for the MHz range**
- **Time resolved optical pulse and power measurements**
- **Laser intensity noise measurements (RIN)**
- **Optical front-end for oscilloscopes, spectrum analyzers, A/D converters and RF lock-in amplifiers**

Block Diagram

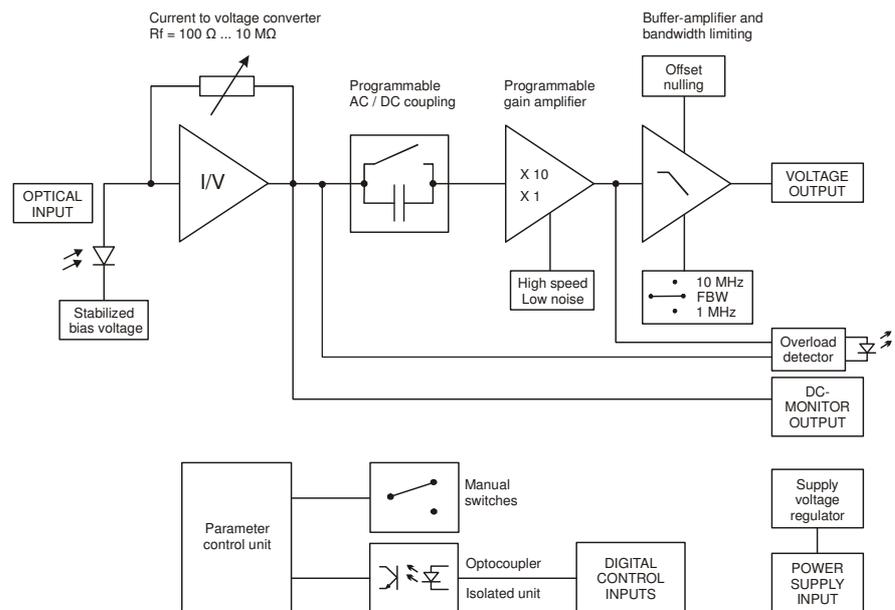

200 MHz Variable Gain Photoreceiver

Available Versions

OE-300-IN-03-FST

1.035"-40 threaded flange
for free space applications and for use with various types
of optical standard accessories

OE-300-IN-03-FS

25 mm dia. unthreaded flange
for free space applications

1.035"-40 threaded flange
Internal threaded coupler ring
with 30 mm outer dia. (included)



OE-300-SI-10-FST

Related OE-300 Models

See separate datasheets for following models on www.femto.de:

OE-300-SI-10-FST

Si-PIN, 1 x 1 mm, 400 - 1000 nm
1.035"-40 threaded flange

OE-300-SI-10-FS

Si-PIN, 1 x 1 mm, 400 - 1000 nm
25 mm dia. unthreaded flange

OE-300-SI-30-FST

Si-PIN, \varnothing 3 mm, 320 - 1000 nm
1.035"-40 threaded flange

OE-300-SI-30-FS

Si-PIN, \varnothing 3 mm, 320 - 1000 nm
25 mm dia. unthreaded flange

OE-300-IN-01-FC

InGaAs-PIN, \varnothing 80 μ m, 900 - 1700 nm
FC fiber receptacle only

OE-300-S

customized versions available on request

Available Accessories

PRA-PAP



post adapter plate,
easy to mount on
FEMTO photoreceiver
series OE, FWPR, HCA-S
and LCA-S

(picture shows model
OE-300-SI-10)



PS-15



power supply,
input: 100 - 240 VAC,
output: \pm 15 VDC, +400/-250 mA

LUCI-10



compact digital I/O interface for USB remote control,
supports opto-isolation of amplifier signal path from PC
USB port, 16 digital outputs, 3 opto-isolated digital inputs,
bus-powered operation

200 MHz Variable Gain Photoreceiver

Specifications

Test conditions $V_s = \pm 15\text{ V}$, $T_A = 25\text{ }^\circ\text{C}$, system impedance = $50\ \Omega$

Gain

Transimpedance gain $1 \times 10^2 \dots 1 \times 10^8\ \text{V/A}$
Gain accuracy $\pm 1\ \%$

Frequency Response

Lower cut-off frequency DC/100 Hz, switchable
Upper cut-off frequency up to 200 MHz (see table below), switchable to 1 MHz or 10 MHz

Input

Noise equivalent power (NEP) see table below
Max. CW saturation power see table below

Detector

Detector InGaAs-PIN photodiode
Active area $300\ \mu\text{m}$ dia.

Spectral response 800 - 1700 nm
Sensitivity R 0.95 A/W typ. @ 1550 nm
Dark current 0.1 nA typ.

Performance Depending on Gain Setting

Gain setting (low noise) (V/A)	10^2	10^3	10^4	10^5	10^6	10^7
Upper cut-off frequency (-3 dB)	200 MHz	80 MHz	14 MHz	3.5 MHz	1.8 MHz	220 kHz
NEP ($\sqrt{\text{Hz}}$, @ 1550 nm)	192 pW	23 pW	1.9 pW	410 fW	152 fW	55 fW
Measured at	20 MHz	8 MHz	1.4 MHz	350 kHz	180 kHz	22 kHz
Integrated input noise (RMS)*	4.8 μW	370 nW	23 nW	3.4 nW	0.82 nW	64 pW
CW sat. power (@ 1550 nm)	10 mW	1.0 mW	100 μW	10 μW	1.0 μW	100 nW
Gain setting (high speed) (V/A)	10^3	10^4	10^5	10^6	10^7	10^8
Upper cut-off frequency (-3 dB)	175 MHz	80 MHz	14 MHz	3.5 MHz	1.8 MHz	220 kHz
NEP ($\sqrt{\text{Hz}}$, @ 1550 nm)	137 pW	6.8 pW	1.4 pW	360 fW	127 fW	52 fW
Measured at	18 MHz	8 MHz	1.4 MHz	350 kHz	175 kHz	22 kHz
Integrated input noise (RMS)*	2.9 μW	270 nW	20 nW	3.3 nW	0.82 nW	64 pW
CW sat. power (@ 1550 nm)	1.0 mW	100 μW	10 μW	1.0 μW	100 nW	10 nW

* The integrated input noise is measured with a shaded input in the full bandwidth ("FBW") setting (referred to 1550 nm). The measurement bandwidth is 3 x the upper cut-off frequency at the specific gain setting; filter slope is a 1st order roll-off.

The input referred peak-peak noise can be calculated from the RMS noise as follows:

$$P_{\text{input noise peak-to-peak}} = P_{\text{input noise RMS}} \times 6$$

The output noise is given by:

$$U_{\text{output noise RMS}} = P_{\text{input noise RMS}} \times \text{gain} \times R$$

$$U_{\text{output noise peak-to-peak}} = U_{\text{output noise RMS}} \times 6 = P_{\text{input noise RMS}} \times \text{gain} \times R \times 6$$

The integrated noise will be reduced considerably by setting the low pass filter to "1 MHz" or "10 MHz" instead of "FBW". This is especially useful for continuous wave (CW) measurements.

200 MHz Variable Gain Photoreceiver

Specifications (continued)

Output	Output voltage range	±1 V (@ 50 Ω load), for linear amplification	
	Output impedance	50 Ω (designed for 50 Ω load)	
Ext. Offset Control	Slew rate	1000 V/μs	
	Max. output current	±40 mA	
	Output offset compensation	adjustable by offset potentiometer and external control voltage, output offset compensation range min. ±100 mV	
	Control voltage range	±10 V	
Indicator LED	Offset control input impedance	15 kΩ	
	Function	overload	
Digital Control	Control input voltage range	LOW bit: -0.8 ... +1.2 V, HIGH bit: +2.3 ... +12 V	
	Control input current	0 mA @ 0 V, 1.5 mA @ +5 V, 4.5 mA @ +12 V	
	Overload output	non active: <0.4 V @ 0 ... -1 mA active: typ. 5 ... 5.1 V @ 0 ... 2 mA	
Power Supply	Supply voltage	±15 V	
	Supply current	+110/-90 mA (depends on operating conditions, recommended power supply capability min ±200 mA)	
	Stabilized power supply output	±12 V, max. 20 mA, +5 V, max. 150 mA	
Case	Weight	320 g (0.74 lb.)	
	Material	AlMg4.5Mn, nickel-plated	
Input Flange	Material	1.4305 stainless steel, glass bead blasted (1.035"-40 threaded flange)	
		AlMg4.5Mn, nickel-plated (25 mm dia. unthreaded flange)	
Coupler Ring	Material	1.4305 stainless steel, glass bead blasted	
DC Monitor Output	Monitor output gain	Mode	Monitor gain
		Low noise	Gain setting divided by -1
	High speed	Gain setting divided by -10	
	Monitor output polarity	inverting	
	Monitor output voltage range	±1 V (@ ≥1 MΩ load)	
	Monitor output bandwidth	DC ... 1 kHz	
	Monitor output impedance	1 kΩ (designed for ≥1 MΩ load)	
Temperature Range	Storage temperature	-40 ... +80 °C	
	Operating temperature	0 ... +60 °C	
Absolute Maximum Ratings	Max. CW power (averaged)	12 mW	
	Digital control input voltage	-5 V/+16 V relative to digital ground DGND (pin 9)	
	Analog control input voltage	±15 V relative to analog ground AGND (pin 3)	
	Power supply voltage	±20 V	

200 MHz Variable Gain Photoreceiver

Remote Control Operation

General

Remote control input bits are opto-isolated and connected by a logical OR function to the local switch settings. For remote control set the corresponding local switches to "Remote", "AC" and "H" and select the desired setting via a bit code at the corresponding digital inputs. Mixed operation, e.g. local AC/DC setting and remote controlled gain setting, is also possible.

Gain setting

Low noise Gain (V/A) Pin 14=HIGH	High speed Gain (V/A) Pin 14=LOW	Pin 12 MSB	Pin 11	Pin 10 LSB
10^2	10^3	LOW	LOW	LOW
10^3	10^4	LOW	LOW	HIGH
10^4	10^5	LOW	HIGH	LOW
10^5	10^6	LOW	HIGH	HIGH
10^6	10^7	HIGH	LOW	LOW
10^7	10^8	HIGH	LOW	HIGH

AC/DC setting

Coupling	Pin 13
DC	LOW
AC	HIGH


Low pass filter setting

Upper cut-off freq. limit	Pin 15	Pin 16
full bandwidth	LOW	LOW
10 MHz	HIGH	LOW
1 MHz	LOW	HIGH

High speed / low noise setting

Mode	Pin 14
low noise mode	LOW
high speed mode	HIGH

Spectral Responsivity

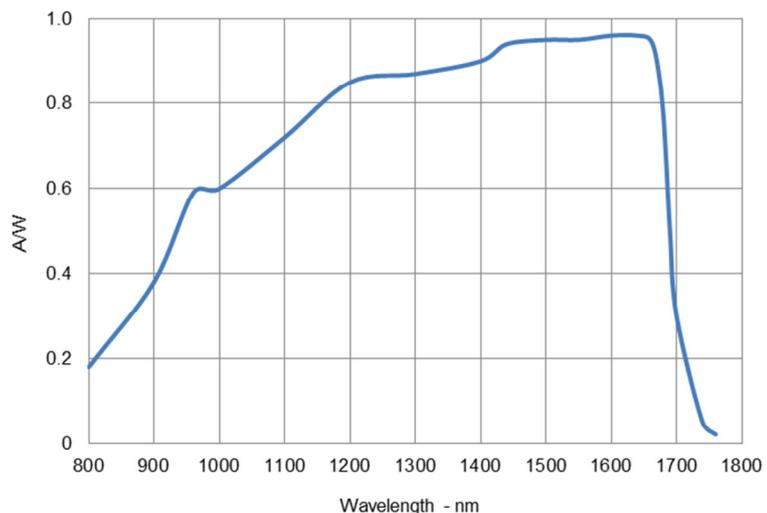


200 MHz Variable Gain Photoreceiver

Typical Performance
Characteristic


Frequency response

$$V_{\text{Supply}} = \pm 15 V_{\text{DC}}; R_{\text{Load}} = 50 \Omega$$


Gain setting: $L10^2, H10^3$

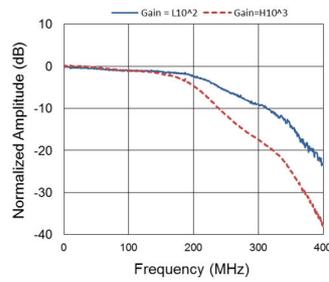


Gain setting: $L10^3, H10^4$

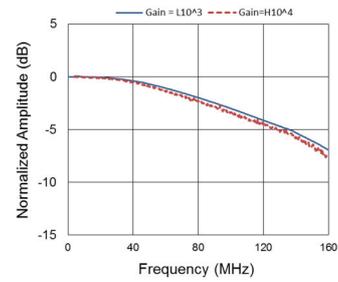


Gain setting: $L10^4, H10^5$

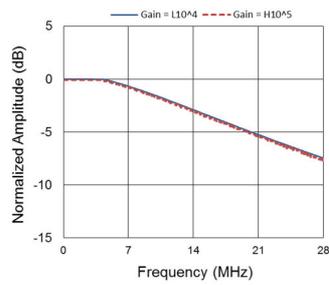


Gain setting: $L10^5, H10^6$

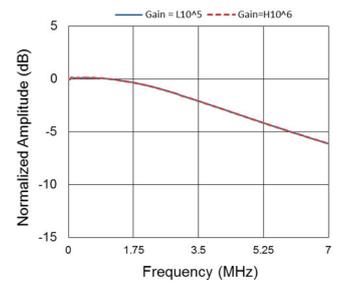

Gain setting: $L10^6, H10^7$

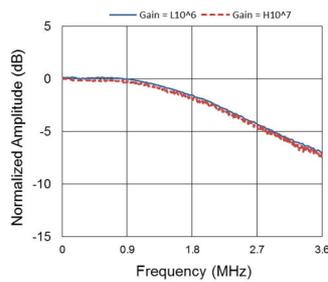

Gain setting: $L10^7, H10^8$

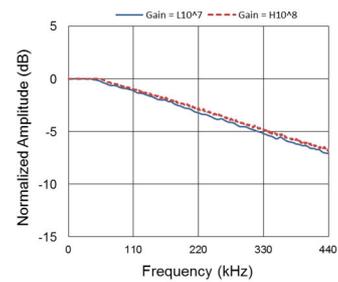


200 MHz Variable Gain Photoreceiver


Typical Performance
Characteristic (continued)

Input noise equivalent power (NEP)


Gain setting $L10^2, H10^3$

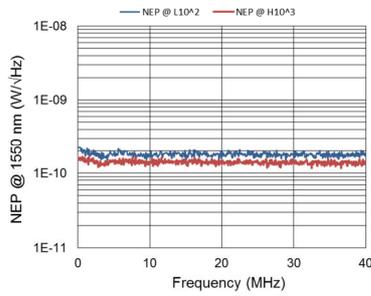


Gain setting $L10^3, H10^4$

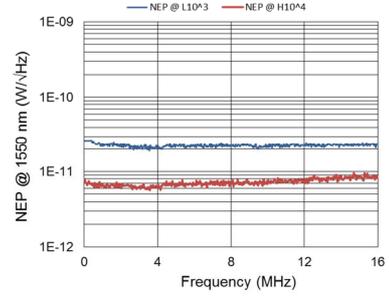


Gain setting: $L10^4, H10^5$

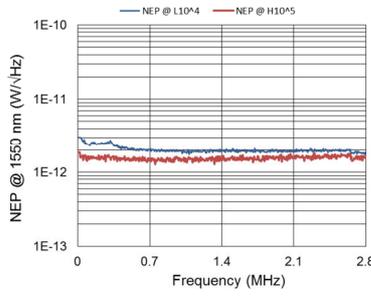


Gain setting: $L10^5, H10^6$

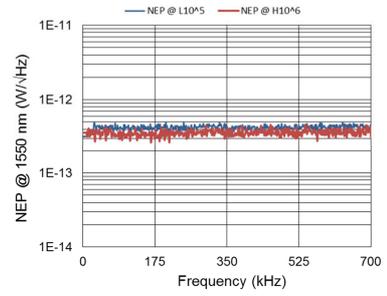

Gain setting: $L10^6, H10^7$

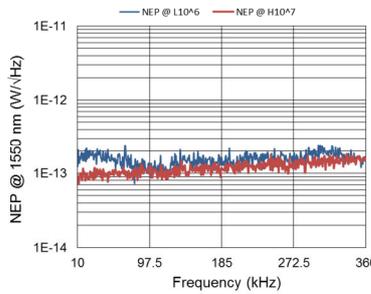

Gain setting: $L10^7, H10^8$

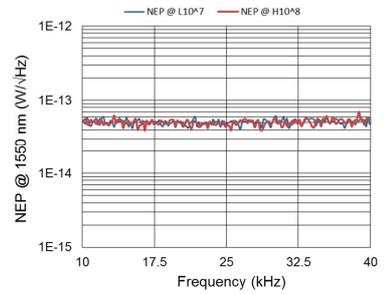


200 MHz Variable Gain Photoreceiver

Typical Performance
Characteristic (continued)


Signal pulse response

Gain setting L10²

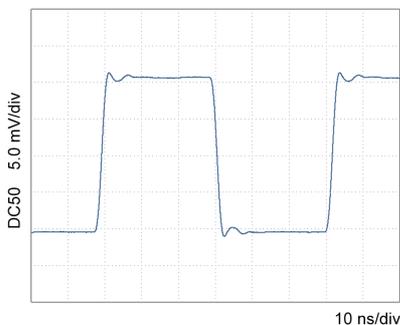


Rise: 1.92 ns Fall: 1.93 ns

Gain setting H10³

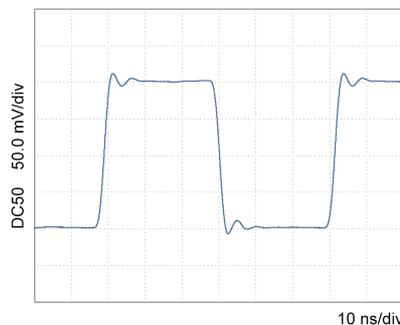


Rise: 2.28 ns Fall: 2.30 ns

Gain setting L10³

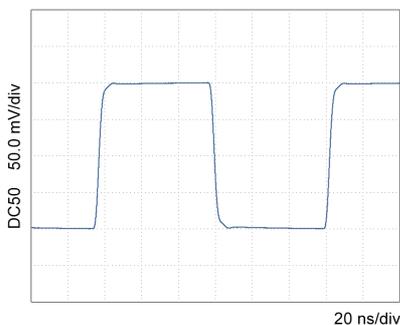


Rise: 3.45 ns Fall: 3.49 ns

Gain setting H10⁴

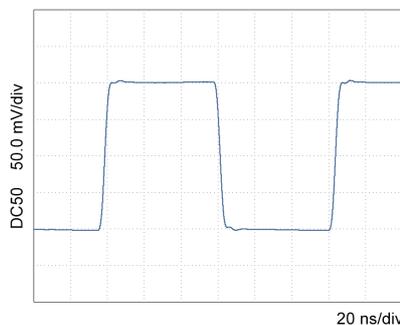


Rise: 3.59 ns Fall: 3.61 ns

Gain setting L10⁴

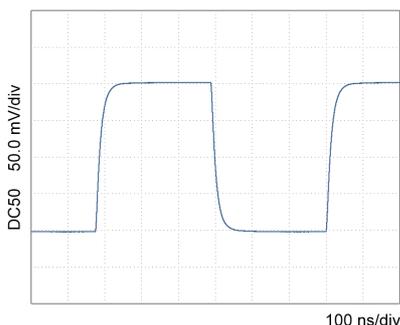

Rise: 25.98 ns Fall: 26.90 ns

Gain setting H10⁵

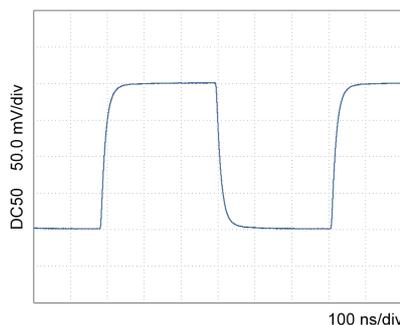


Rise: 27.24 ns Fall: 27.11 ns

200 MHz Variable Gain Photoreceiver


Typical Performance
Characteristic (continued)

Gain setting L10⁵

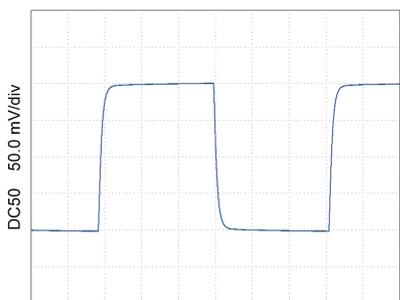


Rise: 82.24 ns Fall: 84.12 ns

Gain setting H10⁶

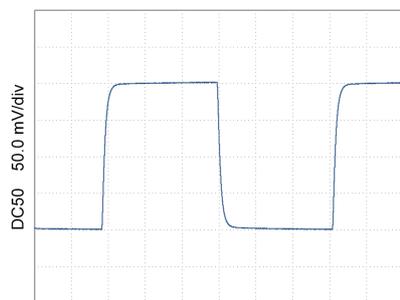

Rise: 81.92 ns Fall: 84.24 ns

Gain setting L10⁶

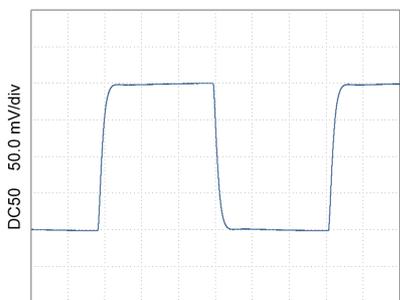


Rise: 208.6 ns Fall: 212.5 ns

Gain setting H10⁷




Rise: 208.8 ns Fall: 208.9 ns

Gain setting L10⁷

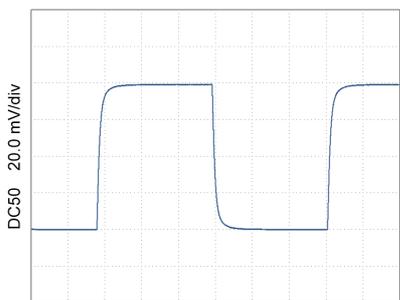

Rise: 1639.2 ns Fall: 1691.2 ns

Gain setting H10⁸

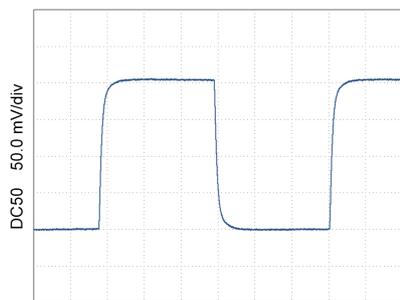


Rise: 1701.6 ns Fall: 1702.4 ns


200 MHz Variable Gain Photoreceiver

Dimensions

Threaded free space input OE-300-IN-03-FST:

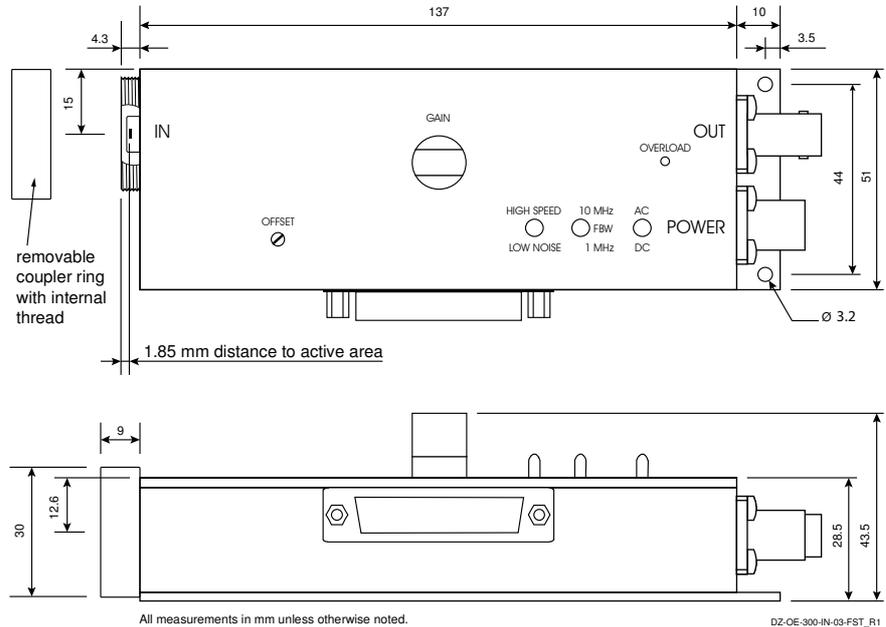

Free space input OE-300-IN-03-FS:

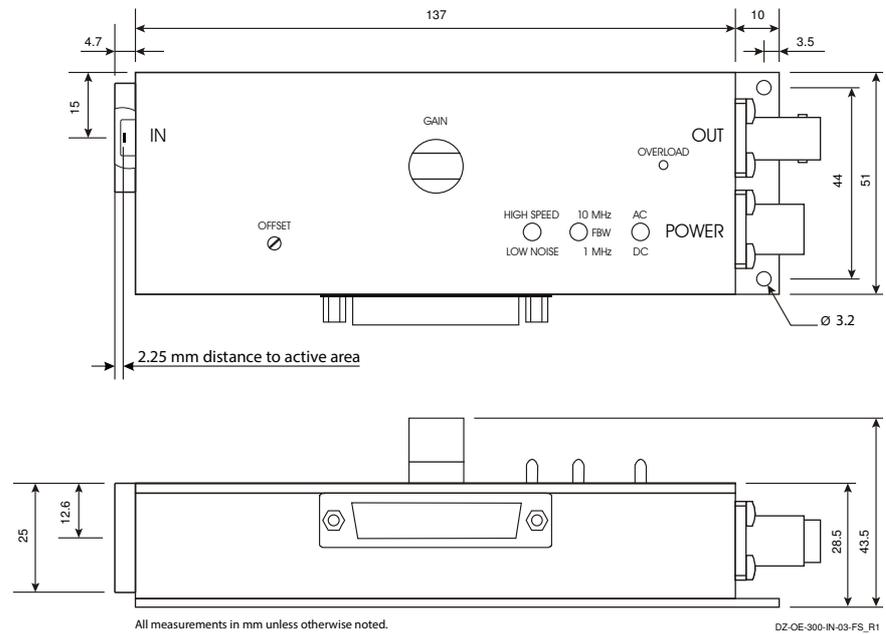

FEMTO Messtechnik GmbH
 Klosterstr. 64
 10179 Berlin · Germany
 Phone: +49 30 280 4711-0
 Fax: +49 30 280 4711-11
 Email: info@femto.de
 www.femto.de

Specifications are subject to change without notice. Information provided herein is believed to be accurate and reliable. However, no responsibility is assumed by FEMTO Messtechnik GmbH for its use, nor for any infringement of patents or other rights of third parties which may result from its use. No license is granted by implication or otherwise under any patent or patent rights of FEMTO Messtechnik GmbH. Product names mentioned may also be trademarks used here for identification purposes only.

© by FEMTO Messtechnik GmbH · Printed in Germany